

## ABgene ALPS300



**Manufacturer:** ABgene  
**Model Number:** Alps300  
**Web Address:** [www.abgene.com](http://www.abgene.com)

The ALPS300 is an automated heat sealer, ideal for medium to high throughput applications. The instrument can seal a wide range of plates, even accommodating different heights.

### Sealer: Commands

- ▶ **S( )** - Seals the microplate.
- ▶ **?( )** - Retrieve the status of the sealer.
- ▶ **C( )** - Retrieves the temperature setpoint (in Celcius).
- ▶ **A( temperature )** - Sets the temperature setpoint (in Celcius) used for sealing.

temperature	Integer	Specifies the temperature (in Celcius) used for sealing. Range of values: <b>100</b> to <b>200</b> .
-------------	---------	--

- ▶ **F( )** - Retrieves the actual temperature (in Celcius) of the sealer.
- ▶ **D( )** - Retrieves the current sealing time in units of 10ths of a second.
- ▶ **B( time )** - Sets the the sealing time in units of 10ths of a second.

time	Integer	Specifies the sealing time. Range of values: <b>0</b> to <b>99</b> .
------	---------	--

- ▶ **E( )** - Retrieve the error code.

### Sealer: Errors

- ▶ **Error( )** - Fatal or non-fatal sealer error.

ReTiSoft Inc.  
 366 Revus Avenue, Unit 21  
 Mississauga, Ontario, Canada, L5G-4S5  
 Main: 647-724-2398 Europe: 33-9-7518-0225  
 Web: [www.retisoft.ca](http://www.retisoft.ca) Email: [prodziew@retisoft.ca](mailto:prodziew@retisoft.ca)