

(c) ReTiSoft Inc.

[www.retisoft.ca](http://www.retisoft.ca)

# Apricot PP150



Manufacturer: Apricot  
 Model Number: Apricot-PP150  
 Web Address: [www.apricotdesigns.com](http://www.apricotdesigns.com)

Apricot PP150 is a unique interchangeable pipetting head system that allows you to use 96, 384 and 1536-well labware on the same machine.

## Dispenser: Commands

▶ **MoveLoadPosition( station )** - Move the loading position.

|         |         |  |
|---------|---------|--|
| station | Integer | Specifies the station to move to. Range of values: 1 to 4. |
|---------|---------|--|

▶ **RunProgram( path )** - Run the program.

|      |        |                                    |
|------|--------|------------------------------------|
| path | String | Specifies the path of the program. |
|------|--------|------------------------------------|

## Dispenser: Errors

🚨 **DispenseError( alarmMessage )** - Error occurred during command execution.

|              |        |                                  |
|--------------|--------|----------------------------------|
| alarmMessage | String | Error message for the instrument |
|--------------|--------|----------------------------------|

ReTiSoft Inc.  
 366 Revus Avenue, Unit 21  
 Mississauga, Ontario, Canada, L5G-4S5  
 Main: 647-724-2398 Europe: 33-9-7518-0225  
 Web: [www.retisoft.ca](http://www.retisoft.ca) Email: [prodziew@retisoft.ca](mailto:prodziew@retisoft.ca)