

## Direct Logic 06 PLC



Manufacturer: Automation Direct  
 Model Number: DL06  
 Web Address: [www.automationdirect.com](http://www.automationdirect.com)

The DL06 micro PLC family has 20 inputs and 16 outputs in the base unit. The DL06 has four option module slots which can be used to add I/O or provide additional communications options.

### DL06 PLC: Commands

▶ **SetOutput( outputLine, state )** - Perform a memory write to the PLC.

outputLine	Integer	Memory location for output from the PLC (1-16). Range of values: 1 to 16.
state	Boolean	Setting the output to true or false.

▶ **GetInput( inputLine )** - Perform a memory read from the PLC (1-20).

inputLine	Integer	Memory location for input from the PLC. Range of values: 1 to 20.
-----------	---------	---

### DL06 PLC: Errors

🔥 **PLCError( description )** - Error occurred during command execution.

description	String	Error description.
-------------	--------	--------------------

ReTiSoft Inc.  
 366 Revus Avenue, Unit 21  
 Mississauga, Ontario, Canada, L5G-4S5  
 Main: 647-724-2398 Europe: 33-9-7518-0225  
 Web: [www.retisoft.ca](http://www.retisoft.ca) Email: [prodziew@retisoft.ca](mailto:prodziew@retisoft.ca)