

Axon ImageXpress 5000A



Manufacturer: Axon Instruments, Inc.
Model Number: ImageXpress-5000A
Web Address: www.axon.com

The ImageXpress 5000A automated cellular imaging and analysis system is an integrated hardware and software system designed specifically for rapid acquisition and analysis of fluorescence images for high throughput cell-based screening.

Cellular Imaging and Analysis System: Commands

▶ **RunScript(script, barcode, plateName)** - Runs the specified script on the instrument.

script	String	Full pathname of the script to run.
barcode	String	Barcode for the current microplate.
plateName	String	Name of the plate for further microplate annotation.

▶ **LoadPlate(position)** - Loads/unloads the microplate into/from the instrument.

position	Integer	Identifies the position for the microplate. 0: microplate is ready to be tended by the robotic arm. 1: microplate is ready for processing.
----------	---------	--

Cellular Imaging and Analysis System: Errors

🔥 **Error(code, description)** - Error occurred during command execution.

code	Integer	Identifies the current error.
description	String	Error description.

ReTiSoft Inc.
 366 Revus Avenue, Unit 21
 Mississauga, Ontario, Canada, L5G-4S5
 Main: 647-724-2398 Europe: 33-9-7518-0225
 Web: www.retisoft.ca Email: prodziew@retisoft.ca