


Beckman Biomek FX

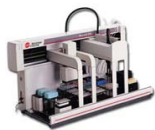

Manufacturer: Beckman Coulter, Inc.
 Model Number: Beckman-Biomek-FX
 Web Address: www.beckmancoulter.com

Biomek FX is a flexible liquid handling system that can simultaneously automate functionality of up to three liquid handlers all on the same worksurface. This liquid handler is offered either preconfigured in smart solutions for targeted applications or custom-tailored to meet your laboratory's individual needs, including the ADMETox Workstation for in vitro permeability analysis of ADME/Tox Assays using the high-throughput method, PAMPA.

Run Methods: Commands

▶ **RunMethod()** - Run the current method on the Biomek FX.

▶ **OpenMethod(methodName)** - Open the specified method on the Biomek FX.

methodName	String	Name of the method to open.
------------	--------	-----------------------------

▶ **GetVariableValue(name)** - Get the value of the specified variable on the Biomek NX.

name	String	Name of the variable.
------	--------	-----------------------

▶ **SetVariableValue(name, value)** - Set the value of the specified variable on the Biomek NX.

name	String	Name of the variable.
------	--------	-----------------------

value	String	Value of the specified variable.
-------	--------	----------------------------------

Pause Methods: Commands

▶ **PauseMethod()** - Pause the current method on the Biomek FX.

▶ **ResumeMethod()** - Resume the current method on the Biomek FX.

Biomek Error

🚫 **BiomekError(code, description)** - Error occurred during command execution.

code	Integer	Error code.
------	---------	-------------

description	String	Error description.
-------------	--------	--------------------

ReTiSoft Inc.
 366 Revus Avenue, Unit 21
 Mississauga, Ontario, Canada, L5G-4S5

Main: 647-724-2398 Europe: 33-9-7518-0225
Web: www.retisoft.ca Email: info@retisoft.ca