

Beckman Biomek NX



Manufacturer: Beckman Coulter, Inc.
 Model Number: Beckman-Biomek-NX
 Web Address: www.beckmancoulter.com

The Biomek NX laboratory automation workstation sets a new standard for flexible laboratory solutions. It puts every aspect of liquid handling - including pipetting, dilution, dispensing, and integrations - into a single, automated system that's as powerful and flexible as it is efficient and economical. The Biomek NX provides all this, while offering unprecedented flexibility for changing needs.

Run Methods: Commands

▶ **RunMethod()** - Run the current method on the Biomek NX.

▶ **OpenMethod(methodName)** - Open the specified method on the Biomek NX.

methodName	String	Name of the method to open.
------------	--------	-----------------------------

▶ **GetVariableValue(name)** - Get the value of the specified variable on the Biomek NX.

name	String	Name of the variable.
------	--------	-----------------------

▶ **SetVariableValue(name, value)** - Set the value of the specified variable on the Biomek NX.

name	String	Name of the variable.
------	--------	-----------------------

value	String	Value of the specified variable.
-------	--------	----------------------------------

Pause Methods: Commands

▶ **PauseMethod()** - Pause the current method on the Biomek NX.

▶ **ResumeMethod()** - Resume the current method on the Biomek NX.

Biomek Error

🔥 **BiomekError(code, description)** - Error occurred during command execution.

code	Integer	Error code.
------	---------	-------------

description	String	Error description.
-------------	--------	--------------------

Web: www.retisoft.ca Email: info@retisoft.ca