

Sagian Plate Carousel



Manufacturer: Beckman Coulter
 Model Number: Beckman-Carousel
 Web Address: www.beckmancoulter.com

SAGIAN Hotels provide plate and tip storage and ambient incubation for your automated assay system. When large unattended capacity is desired, a hotel carousel is a space-saving option. It allows true random access for up to ten hotels containing as many as 18 standard microplates. The carousel rotates, enabling any one of the ten hotels in the structure to be accessed by the robotic arm.

Carousel: Commands

▶ **Home()** - Home the carousel and move to the hotel position defined as position 1.

▶ **Advance(hotels)** - Move the carousel forward the specified number of hotels.

hotels	Integer	Specifies the number of hotels to move forward.
--------	---------	-------------------------------------------------

▶ **MoveToHotel(hotel)** - Move the carousel to the specified hotel.

hotel	Integer	Specifies the hotel number.
-------	---------	-----------------------------

Carousel: Errors

🔥 **CarouselError(alarmMessage)** - Error occurred during command execution.

alarmMessage	String	Error message for the instrument
--------------	--------	----------------------------------

ReTiSoft Inc.
 366 Revus Avenue, Unit 21
 Mississauga, Ontario, Canada, L5G-4S5
 Main: 647-724-2398 Europe: 33-9-7518-0225
 Web: www.retisoft.ca Email: info@retisoft.ca