

Beckman CS Incubator



Manufacturer: Beckman Coulter, Inc.
Model Number: Beckman-CS-Incubator
Web Address: www.beckmancoulter.com

The CS Incubator (4- or 8-position) provides microplate storing capability at temperatures between 4 - 45C using an appropriate circulating temperature bath. The 8-position incubator has a capacity of eight lidded or unlidded standard microplates. The 4-position incubator has a capacity of four lidded or unlidded standard microplates or deep-well microplates or reservoirs.

CS Incubator: Commands

▶ **GetStatus()** - Retrieve the current status of the incubator.

▶ **SetApproachTime(time)** - Set the length of time, in milliseconds, a tray moves at approach speed.

| | | |
|------|---------|---------------------------------|
| time | Integer | Length of time in milliseconds. |
|------|---------|---------------------------------|

▶ **SetApproachSpeed(speed)** - Set the approach speed of the trays.

| | | |
|-------|---------|------------------------------|
| speed | Integer | Approach speed of the trays. |
|-------|---------|------------------------------|

▶ **SetRunTime(time)** - Set the length of time, in milliseconds, a tray moves at run speed.

| | | |
|------|---------|---------------------------------|
| time | Integer | Length of time in milliseconds. |
|------|---------|---------------------------------|

▶ **SetRunSpeed(speed)** - Set the run speed of the trays.

| | | |
|-------|---------|-------------------------|
| speed | Integer | Run speed of the trays. |
|-------|---------|-------------------------|

▶ **MoveTray(tray, move)** - Move a tray in or out of the incubator.

| | | |
|------|---------|--|
| tray | Integer | Tray to be moved. Range of values: 1 to 8. |
|------|---------|--|

| | | |
|------|---------|---|
| move | Boolean | True to move the tray in, false to move the tray out. |
|------|---------|---|

▶ **ResetTrays()** - Moves all the trays into the incubator.

CS Incubator: Errors

🔥 **IncubatorError(code)** - Error occurred during command execution.

| | | |
|------|--------|-------------|
| code | String | Error code. |
|------|--------|-------------|