

(c) ReTiSoft Inc.

www.retisoft.ca

Beckman Forma Incubator



Manufacturer: Beckman Coulter, Inc.
 Model Number: Beckman-Forma-Incubator
 Web Address: www.beckmancoulter.com

Forma Scientific Environmental CO2 Incubator features bright and clear digital displays, temperature control dials and a rotating carousel.

Forma Incubator: Commands

▶ **Home()** - Home the incubator and move to the hotel position defined as position 1.

▶ **Advance(hotels)** - Move the incubator forward the specified number of hotels.

hotels	Integer	Specifies the number of hotels to move forward.
--------	---------	---

▶ **MoveToHotel(hotel)** - Move the incubator to the specified hotel.

hotel	Integer	Specifies the hotel number.
-------	---------	-----------------------------

▶ **DoorOpen()** - Open the incubator door.

▶ **DoorClose()** - Close the incubator door.

Incubator Error

🔥 **IncubatorError(code)** - Error occurred during command execution.

code	String	Error code.
------	--------	-------------

ReTiSoft Inc.
 366 Revus Avenue, Unit 21
 Mississauga, Ontario, Canada, L5G-4S5
 Main: 647-724-2398 Europe: 33-9-7518-0225
 Web: www.retisoft.ca Email: info@retisoft.ca