


## Brandel RS3000

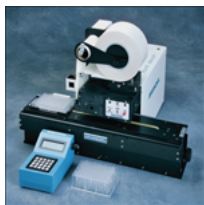

**Manufacturer:** Brandel Biosciences  
**Model Number:** RS3000  
**Web Address:** [www.brandel.com](http://www.brandel.com)

Brandel's plate sealer offers a pressure-based solution for applying a uniform siliconized or adhesive seal for most any plate configuration, height or composition. The sealer can be used with a variety of sealing tapes, including transparent, translucent, foils and gas permeable. A variety of sealing tapes are suitable for use with most chemicals (including DMSO), compatible with cells, and capable of withstanding temperatures as low as -80C.

### Plate Sealer: Commands

▶ **XMLP( profile )** - Load and seal a plate with the profile stored in memory. Return the plate to the home position.

profile	Integer	Number of the profile stored in memory. Range of values: <b>0</b> to <b>9</b> .
---------	---------	---

▶ **XMLp( profile )** - Load and seal a plate with the profile stored in memory. Put the plate on the other side.

profile	Integer	Number of the profile stored in memory. Range of values: <b>0</b> to <b>9</b> .
---------	---------	---

### Plate Sealer: Errors

🔥 **SealerError( message )** - Error occurred on the sealer.

message	String	Error message.
---------	--------	----------------

ReTiSoft Inc.  
 366 Revus Avenue, Unit 21  
 Mississauga, Ontario, Canada, L5G-4S5  
 Main: 647-724-2398 Europe: 33-9-7518-0225  
 Web: [www.retisoft.ca](http://www.retisoft.ca) Email: [prodziew@retisoft.ca](mailto:prodziew@retisoft.ca)