

Computype LAP4110



Manufacturer: Computype
Model Number: LAP4110
Web Address: www.computype.com

The LAP4110 is a versatile print and apply label applicator that is built on an industrial Zebra thermal transfer label printer. It features a zero-queue, fixed position horizontal swing tamp applicator which places labels with extreme accuracy (+/- .020 in.) and reliability.

Run - Label Applicator: Commands

▶ **OpenJob(path)** - Opens a job without displaying the browse dialog box.

path	String	Full file path of an existing and valid job.
------	--------	--

▶ **CloseJob()** - Closes the job record that is currently being referenced.

▶ **Start()** - Prepares job record and applicator for printing, applying and scanning operation. Must be executed before the PrintJob command.

▶ **Stop()** - Closes applicator for printing, applying and scanning operation.

▶ **UseExternalSerialNumber(flag)** - Determines whether to use a serial number generated by the job record, or a value supplied by the client application for printing labels.

flag	Boolean	true: use a serial number supplied by the client application, false: use internal numbers generated by the job record.
------	---------	---

▶ **SetSerialNumber(serialNumber)** - Sets the string value of the current serial number.

(Note: If UseExternalSerialNumber is true then serial number value is expected to be supplied by the client application).

serialNumber	String	Serial number to be printed.
--------------	--------	------------------------------

▶ **PrintJob()** - Prints, applies and optionally scans a label based on the name of the job file currently open.

▶ **Status()** - Retrieves an integer value that corresponds to the instrument status. 0: A general type error has occurred. 1: Job record is loaded, applicator has been initialized. 2: Wait for a part to be detected. 4: Label template has been sent to printer, waiting for the cycle complete signal from the Morphor3. 8: Label has been applied, waiting for the part to be removed. 16: Label has been printed, applied and scanned successfully. 32: Label has not been scanned because of the software configuration. 64: Barcode scanned is not the same as the barcode printed. 128: Barcode was scanned successfully. 256: The job has stopped, the serial communication to the Morphor3 has been closed, and the applicator is at its rest state. 521: Error connecting to the Morphor3. 1024: Applicator is not ready to access printer templates. 2048: Error applying a label. 5000: Waiting for scan operation to complete. 5001: Error occurred during the scan operation. 5002: Scan has completed successfully. 6000: Creating a new job thread in the JoTS application server.

Run - Label Applicator: Errors

🔥 **Error(number)** - Fatal or non-fatal LAP4110 error has occurred.

number	Integer	Number of the error.
--------	---------	----------------------

ReTiSoft Inc.
 366 Revus Avenue, Unit 21
 Mississauga, Ontario, Canada, L5G-4S5
 Main: 647-724-2398 Europe: 33-9-7518-0225
 Web: www.retisoft.ca Email: prodziew@retisoft.ca