


Dynamic Devices Nectar

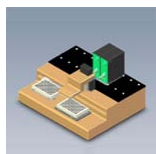

Manufacturer: Dynamic Devices
 Model Number: Dynamic-Devices-Nectar
 Web Address: www.dynamicdevices.com

Nectar, the Independent Channel Bulk Reagent Dispenser, handles 96 or 384 well plates and can be configured with up to 8 independent reagent modules. Any single reagent can be dispensed at any given time.

Nectar Dispenser: Commands

▶ **ShowWindow()** - Show the Method Manager window.

▶ **HideWindow()** - Hide the Method Manager window.

▶ **StartMethod(methodName)** - Start the specified method on the Nectar.

methodName	String	Path for the specified method.
------------	--------	--------------------------------

▶ **CloseMethod(methodName)** - Close the specified method on the Nectar.

methodName	String	Path for the specified method.
------------	--------	--------------------------------

▶ **SetVariable(variableName, variableValue)** - Sets the value for specified variable.

variableName	String	Name of the variable.
--------------	--------	-----------------------

variableValue	String	Value for the variable.
---------------	--------	-------------------------

Nectar Dispenser: Errors

🔥 **OasisError(code, description)** - Error occurred during command execution.

code	Integer	Error code.
------	---------	-------------

description	String	Error description.
-------------	--------	--------------------

ReTiSoft Inc.
 366 Revus Avenue, Unit 21
 Mississauga, Ontario, Canada, L5G-4S5
 Main: 647-724-2398 Europe: 33-9-7518-0225
 Web: www.retisoft.ca Email: prodziew@retisoft.ca