

## Hamilton Microlab4200-96MPH



**Manufacturer:** Hamilton Company  
**Model Number:** Microlab4200-MPH96  
**Web Address:** [www.hamiltoncompany.com](http://www.hamiltoncompany.com)

The Microlab4200 MPH-96 is a high throughput liquid handling workstation with 96 pipetting channels and syringes for optimal performance and accuracy. It also contains high efficiency fixed probe wash system.

### Liquid Handler: Commands

▶ **RunMethod( )** - Run the currently set method on the instrument.

▶ **SetDeckLayoutName( layoutFile )** - Set the name of the deck layout file.

layoutFile	String	name of the deck layout file.
------------	--------	-------------------------------

▶ **SetMethodName( methodName )** - Set the name of the method file.

methodName	String	name of the method to be executed on the instrument.
------------	--------	--

▶ **SetTraceName( traceFile )** - Set the name of the trace file.

traceFile	String	name of the trace file.
-----------	--------	-------------------------

▶ **Enable( enable )** - Enable/disable the remote control of the instrument.

enable	Boolean	enable/disable flag for remote control of the instrument.
--------	---------	---

### Liquid Handler: Errors

🔥 **MethodTimeout( )** - Method could not be invoked. Timeout period has expired.

🔥 **MethodAborted( )** - Method was aborted.

🔥 **MethodFailed( code )** - An error occurred during method execution.

code	Integer	Error code.
------	---------	-------------

ReTiSoft Inc.  
 366 Revus Avenue, Unit 21  
 Mississauga, Ontario, Canada, L5G-4S5

Main: 647-724-2398 Europe: 33-9-7518-0225

Web: [www.retisoft.ca](http://www.retisoft.ca) Email: [prodziew@retisoft.ca](mailto:prodziew@retisoft.ca)