

Myomics M.A.D.



Manufacturer: Myomics Inc.
 Model Number: Myomics-MAD
 Web Address: www.myomics.com

The Myoforce Assay Device (M.A.D.) is able to electrically stimulate and capture force data from muscle tissues.

Run: Commands

- ▶ **EjectPlate()** - Open the door and eject the plate.
- ▶ **AcceptPlate()** - Close the door and accept the plate.
- ▶ **RunMeasurement(measurementID)** - Run the specified measurement.

measurementID	String	Measurement identifier.
---------------	--------	-------------------------

- ▶ **Clean(slotID)** - Clean the plate positioned on the specified slot.

slotID	Integer	Slot identifier. Range of values: 1 to 10.
--------	---------	--

- ▶ **EjectClean(slotID)** - Open the door and remove the plate from the specified slot.

slotID	Integer	Slot identifier. Range of values: 1 to 10.
--------	---------	--

- ▶ **AcceptClean(slotID)** - Close the door and move the plate onto the specified slot.

slotID	Integer	Slot identifier. Range of values: 1 to 10.
--------	---------	--

Run: Errors

- 🔥 **ReaderError(alarmCode)** - Error occurred during command execution.

alarmCode	Integer	Error code for the reader.
-----------	---------	----------------------------

Abort: Commands

- ▶ **Shutdown()** - Stop the measurement, eject the plate and shutdown the instrument.

Abort: Errors

- 🔥 **ReaderError(alarmCode)** - Error occurred during command execution.

alarmCode	Integer	Error code for the reader.
-----------	---------	----------------------------

ReTiSoft Inc.
366 Revus Avenue, Unit 21
Mississauga, Ontario, Canada, L5G-4S5
Main: 647-724-2398 Europe: 33-9-7518-0225
Web: www.retisoft.ca Email: prodziew@retisoft.ca