

PerkinElmer PlateStak



Manufacturer: PerkinElmer Life Sciences
 Model Number: PerkinElmer-PlateStak
 Web Address: www.perkinelmer.com

The PerkinElmer® PlateStak is a superior solution for addressing increased capacity and unattended processing needs by automating handling and storage of microplates and RoboRack tip boxes.

PlateStak: Commands

- ▶ **ScanBarcode()** - Turn on the barcode scanner and scan a barcode.
- ▶ **GetLocation()** - Get the current location of the shuttle.
- ▶ **MoveTray(source, destination)** - Move one tray from source to destination.

source	Integer	Specifies the source location. Range of values: 2 to 3.
destination	Integer	Specifies the destination location. Range of values: 1 to 4.

- ▶ **MoveTrayWithScan(source, destination)** - Move one tray from source to destination with barcode scanning.

source	Integer	Specifies the source location. Range of values: 2 to 3.
destination	Integer	Specifies the destination location. Range of values: 1 to 4.

PlateStak: Errors

- 🔥 **HandlerError(alarmMessage)** - Error occurred during command execution.

alarmMessage	String	Error message for the instrument
--------------	--------	----------------------------------

ReTiSoft Inc.
 366 Revus Avenue, Unit 21
 Mississauga, Ontario, Canada, L5G-4S5
 Main: 647-724-2398 Europe: 33-9-7518-0225
 Web: www.retisoft.ca Email: info@retisoft.ca