

Tecan Columbus



Manufacturer: Tecan
Model Number: Columbus
Web Address: www.tecan.com

Columbus washer is a flexible, fully automated microplate strip washer for 96- and 384-well plates.

Plate Washer: Commands

- ▶ **Y0()** - Transfer the washer into the remote control.
- ▶ **E()** - Transfer the washer back into the manual (local) control.
- ▶ **S(pl, p2)** - Start the program on the washer.

pl	Integer	Number of the program to start.
p2	Integer	Number of the strips on the plate.

Plate Washer: Errors

🔥 **WasherError(alarmCode)** - Error occurred during command execution.

alarmCode	Integer	Error code: 1 - init error, 2 - invalid command, 3 - invalid operand, 4 - invalid command sequence, 5 - system error (not used), 6 - res, 7 - plate mismatch, 8 - plate transport error, 9 - manifold transport error, 10 - dispenses pump error, 11 - fill verification error, 12 - manifold error, 13 - waste full error, 14 - test not defined, 15 - command overflow, 16 - fill verification function error, 17 - fill verification aspiration error.
-----------	---------	---

ReTiSoft Inc.
 366 Revus Avenue, Unit 21
 Mississauga, Ontario, Canada, L5G-4S5
 Main: 647-724-2398 Europe: 33-9-7518-0225
 Web: www.retisoft.ca Email: prodziew@retisoft.ca