

## Tecan Sunrise



**Manufacturer:** Tecan  
**Model Number:** Sunrise  
**Web Address:** [www.tecan.com](http://www.tecan.com)

Sunrise combines fast, reproducible and accurate measurements with high-quality optical performance for diagnostic and research laboratories. Sunrise supports a large variety of absorbance based immunoassay protocols (immunometric methods, such as ELISA assays; competitive and homo- or heterogenous immunoassays), kinetic assays or colorimetric cell toxicity assays amongst numerous further applications.

### Plate Reader: Commands

▶ **GetServerVersion( )** - Retrieve the version and licence information for the Magellan server.

▶ **PlateIn( )** - Move the plate carrier into the instrument.

▶ **PlateOut( )** - Move the plate carrier out of the instrument.

▶ **GetPlatePosition( )** - Retrieve the current plate position.

▶ **SetFileName( fileName )** - Set the name of the output file for the method.

fileName	String	Name of the output file for the method.
----------	--------	---

▶ **SetEmpty( state )** - Reset the state of the workspace.

state	Boolean	True resets the workcell to the empty state.
-------	---------	--

▶ **InsertMethod( method )** - Load the specified method into the current workspace.

method	String	Name of the method to be loaded.
--------	--------	----------------------------------

▶ **Run( )** - Run the previously loaded method.

### Plate Reader: Errors

🔥 **SunriseError( description )** - Error occurred during command execution.

description	String	Error description
-------------	--------	-------------------

ReTiSoft Inc.  
366 Revus Avenue, Unit 21  
Mississauga, Ontario, Canada, L5G-4S5  
Main: 647-724-2398 Europe: 33-9-7518-0225  
Web: [www.retisoft.ca](http://www.retisoft.ca) Email: [prodziew@retisoft.ca](mailto:prodziew@retisoft.ca)