

## Tecan ULTRA Evolution



**Manufacturer:** Tecan  
**Model Number:** ULTRA Evolution  
**Web Address:** [www.tecan.com](http://www.tecan.com)

The Ultra Evolution is a multifunctional high performance detection platform. Following the need for having access to most applications in one easy-to-use instrument the Ultra Evolution offers to the user up to nine detection modes in one compact instrument. Flexibility, high speed and sensitivity as well as reliability open a wide field of application areas including high throughput screening, assay development, secondary and confirmatory screening.

### Plate Reader: Commands

- ▶ **GetServerVersion( )** - Retrieve the version and licence information for the Magellan server.
- ▶ **PlateIn( )** - Move the plate carrier into the instrument.
- ▶ **PlateOut( )** - Move the plate carrier out of the instrument.
- ▶ **GetPlatePosition( )** - Retrieve the current plate position.
- ▶ **SetFileName( fileName )** - Set the name of the output file for the method.

fileName	String	Name of the output file for the method.
----------	--------	---

- ▶ **SetEmpty( state )** - Reset the state of the workspace.

state	Boolean	True resets the workcell to the empty state.
-------	---------	--

- ▶ **InsertMethod( method )** - Load the specified method into the current workspace.

method	String	Name of the method to be loaded.
--------	--------	----------------------------------

- ▶ **Run( )** - Run the previously loaded method.

### Plate Reader: Errors

- 🔥 **ULTRA EvolutionError( description )** - Error occurred during command execution.

description	String	Error description
-------------	--------	-------------------

ReTiSoft Inc.  
366 Revus Avenue, Unit 21  
Mississauga, Ontario, Canada, L5G-4S5  
Main: 647-724-2398 Europe: 33-9-7518-0225  
Web: [www.retisoft.ca](http://www.retisoft.ca) Email: [prodziew@retisoft.ca](mailto:prodziew@retisoft.ca)