

Tomtec Harvester9600



Manufacturer: Tomtec Inc.
Model Number: Harvester9600
Web Address: www.tomtec.com

Harvester9600 contains removable wash tips, "hot" and "cold" waste, and can be controlled remotely or work as a stand alone unit. The Harvester 9600 is available in filter plate or filter mat formats for robotic operation. Standard glass fiber filter mats for the Betaplate or MicroBeta are carried in frames for robotic handling. UniFilter plates for the Topcount or MicroBeta are handled directly. Special Hotels are available for either format.

Plate Washer: Commands

▶ **gv()** - Retrieves the model and software version numbers.

▶ **ru(N)** - Runs the specified user program.

N	Integer	Specifies the user program. Range of values: 1 to 40 .
---	---------	--

▶ **rp()** - Resumes the execution of the program until the program ends or it is paused.

▶ **st()** - Stops the execution of the previously paused program.

Plate Washer: Errors

🔥 **error(alarmCode)** - Error occurred during command execution.

alarmCode	String	Error code which describes the type of instrument malfunction.
-----------	--------	--

ReTiSoft Inc.
 366 Revus Avenue, Unit 21
 Mississauga, Ontario, Canada, L5G-4S5
 Main: 647-724-2398 Europe: 33-9-7518-0225
 Web: www.retisoft.ca Email: prodziew@retisoft.ca