

Tomtec Quadra-Dispenser



Manufacturer: Tomtec Inc.
Model Number: Quadra-Dispenser
Web Address: www.tomtec.com

The Quadra 384 is a true 384-well pipettor aspirating or dispensing to 384 wells simultaneously. Each piston moves within its' own separate tip in a positive displacement mode, permitting precision pipetting at sub microliter volumes. The three position shuttle is ideal for low volume protocols or assay development.

Liquid Handler: Commands

▶ **qv()** - Retrieves the model and software version numbers.

▶ **gb()** - Returns data from the bar code reader.

▶ **ru(N)** - Runs the specified user program.

N	Integer	Specifies the user program. Range of values: 1 to 40 .
---	---------	--

▶ **rp()** - Resumes the execution of the program until the program ends or it is paused.

▶ **st()** - Stops the execution of the previously paused program.

Liquid Handler: Errors

🔥 **error(alarmCode)** - Error occurred during command execution.

alarmCode	String	Error code which describes the type of instrument malfunction.
-----------	--------	--

ReTiSoft Inc.
 366 Revus Avenue, Unit 21
 Mississauga, Ontario, Canada, L5G-4S5
 Main: 647-724-2398 Europe: 33-9-7518-0225
 Web: www.retisoft.ca Email: prodziew@retisoft.ca